

USER MANUAL

# LED Gooseneck Illuminator

For all AMPS-ILED-23



## Setup

1. Remove illuminator and AC power supply from packaging.
2. Place illuminator on flat, solid surface.

## Operation

1. Connect the AC power adapter to the input jack located on the back of the illuminator base.
2. Plug AC power adapter into wall socket.
3. Switch the main power switch to the ON "I" position.
4. Place the "CHARGE/ON/OFF" Lever in the desired position. (Left = off and charging battery / Middle = On / Right = off and not charging battery)
5. Turn the Illumination intensity knob to the desired intensity
6. Adjust goosenecks to desired position to illuminate sample.



**LAXCO Inc.**

18303 Bothell-Everett Hwy , Suite 140  
Mill Creek, Washington 98012

TEL: +1 425-686-3081

EMAIL: [info@laxcoinc.com](mailto:info@laxcoinc.com)

## SPECIFICATIONS

Dimensions	9.0cm x 10.8cm x 38.0cm
Weight	1.10kg
Operating and Storage Temperature Range	-30°C to 85°C
Relative Humidity	70% or less
Lamp Type	Ultra-bright LED
LED Lifetime	50,000 hours
Color Temperature	approx. 6,800°K
Spectrum Range	White (450nm-750nm)
Power	White: 3W
Power Supply	AC Adapter; Input 100-240V, 50-60Hz Output 9V, 0.6A max

**Statement of Limited Product Warranty:** This model is warranted to be free of all defects in material and workmanship for a period of 12 months (1 year) from the date of delivery. The warranty does not apply to any instrument which has become worn, defective, damaged or broken due to abuse, misuse, tampering, or unauthorized repairs. Under this warranty, Laxco will repair or replace, without charge to the purchaser, any part which upon our examination, appears to be defective in materials or workmanship. Returned Goods Policy for Repair or Replacement Parts: To return goods for repair or replacement, please contact Laxco Customer Service.

**Disclaimer:** Information provided by Laxco is believed to be accurate and reliable. However, no responsibility is assumed by Laxco for its use. The information contained in this publication is subject to change without notice. Words that Laxco considers trademarks have been identified as such. However, neither the presence nor absence of such identification affects the legal status of any trademark. Units of measurement in this document conform to SI standards and practices. All rights reserved. No part of this publication may be reproduced, stored in a retrieval system, or transmitted in any form, be it electronically, mechanically, or by any other means such as photocopying, recording, or otherwise, without the prior written permission of Laxco Inc.



**LAXCO Inc.**

18303 Bothell-Everett Hwy , Suite 140  
Mill Creek, Washington 98012

TEL: +1 425-686-3081 | EMAIL: [info@laxcoinc.com](mailto:info@laxcoinc.com)